


QUICK START GUIDE: SERIES 12000 BALANCES


1. Unpack Balances

- Do not drop balance or orient upside down or on its sides
- Do not push sideways on installed weighing pan

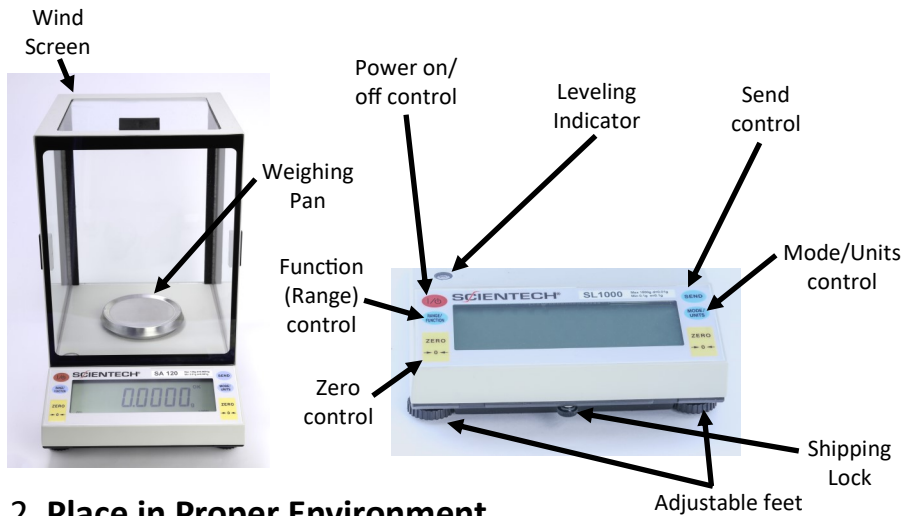


2. Place in Proper Environment

- Operate indoors in a dry environment free from air flow, vibration sources, variations in temperature (no greater than $\pm 10^{\circ}\text{C}$), and with no static electricity.

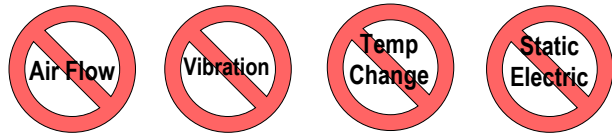


3. Level Balance

- Adjust front feet to center the leveling bubble indicator

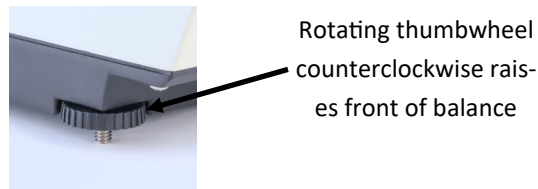

4. Release Shipping Lock


- **Must release on all models for balance to work**
- **Install weighing pan**

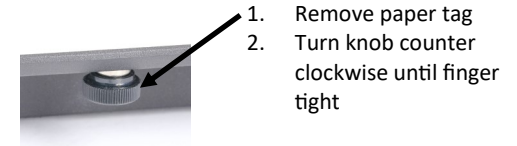

5. Connect Balance to Electric Power

- Make sure AC power adapter voltage and plug type are compatible with local AC power
- Ensure power cord is not a tripping hazard
- Disconnect balance from power when cleaning

6. Power On Balance

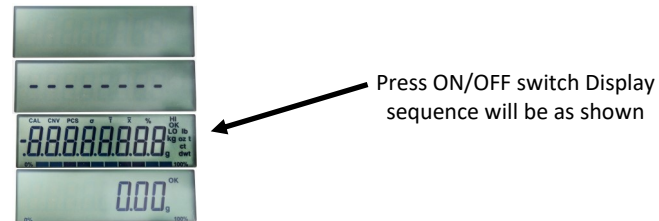

7. Allow 1 hour Warmup and Then Calibrate

- Calibration for internal weight balance: Refer to Operating Manual
- Calibration using external weight: Refer to Operating Manual

8. Where to Get Operating Manual

Download at: www.scientech-inc.com/event-knowledge-center/knowledge-center/quick-start-guides.

9. Support

- If operating support is needed call Scientech Inc. at:
Tel: 303 444 1361 or 1 800 525 0522 or
Email: inst@scientech-inc.com
- If repair is required, contact authorized Scientech repair center. Disassembling balance by non-authorized repair center voids warranty.